USB-Mikroskop-Kamera 500fach

Best.-Nr. 19 13 89

Version 11/11

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Item no. 19 13 89

Version 11/11
CE

Use in accordance with manufacturer's instructions

The product is a camera designed to transfer the recorded image via a USB port to a PC. Thanks to the integrated lens, particularly small objects/surfaces can be magnified and imaged. The power is supplied via a USB port.

This product complies with the applicable National and European requirements. All names of companies and products are the trademarks of the respective owners. All rights reserved.

Package Contents

- Microscope camera
- Table stand
- CD with driver/software
- Operating Instructions

Safety Instructions

The warranty will be void in the event of damage caused by failure to observe these safety instructions! We do not assume any liability for any consequential damage!

We do not assume any liability for material and personal damage caused by improper use or non-compliance with the safety instructions. In such cases the warranty will be null and void.

- The unauthorised conversion and/or modification of the product is not permissible for reasons of safety and regulation (CE).
- The product should only be used in dry indoor areas, it must not be exposed to damp or wet!
- The product is not a toy and should be kept out of the reach of children.
- **Caution, LED light:**
Do not look directly into the LED light!
Do not look into the beam directly or with optical instruments!
- Do not leave packing materials unattended. They may become dangerous playthings for children.
- Handle the product with care; it can be damaged by impacts, blows, or accidental drops, even from a low height.

Driver/Software Installation

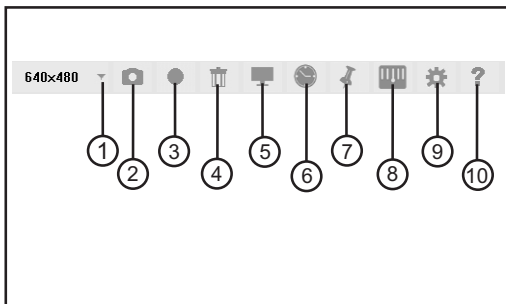
Operating system and application software are subject to permanent change by updates. Therefore not all contingencies can be pointed out in these operating instructions, but solely the basic software installation.

- Start your operating system. Place the CD included in the delivery in an appropriate drive on your computer.
- The software now starts automatically and you can install the software. If the CD does not start automatically, you must start the installation program on the CD manually (main directory on the CD, e.g. "Autorun.exe" etc.). Follow all of the software's instructions. The serial code required by a dialogue window can be found on the CD cover.

On the desktop of your computer a new icon is created. Now connect the microscopic camera to a free USB2.0 port on your computer. The operating system recognises the new hardware and completes the driver installation. Please read the manufacturer's information that can be found on the CD.


- Start the software by double clicking on the newly created icon.
- A new window is opened. Now the program and the micro camera are ready for operation.
- To operate the software different symbols will be shown:

- | | |
|-----------------------|--------------------|
| 1 Resolution settings | 2 Create photo |
| 3 Video recording | 4 Delete files |
| 5 Full screen | 6 Date and time |
| 7 "Freeze" image | 8 Measurement grid |
| 9 Settings | 10 "Help file" |

Controls

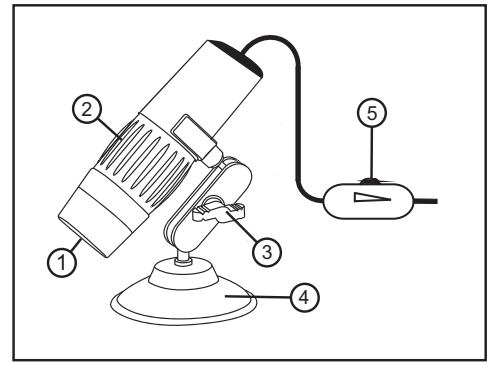- 1 Objective (with LED illumination)
- 2 Adjustment dial
- 3 Locking bolt
- 4 Stand
- 5 Rotary knob (LED brightness)

Operation

- By means of the rotary knob in the connection cable (5) the LED illumination in the front of the lens can be turned off or the brightness can be adjusted steplessly.
- Move the lens (1) of the microscope camera as close as possible to the desired object (e. g. place the front on a newspaper).
- Adjust the picture until it the image is sharp by using the adjustment dial (2).

Do not overwind the adjustment dial. Turn the adjustment wheel slowly, so that you can focus the picture accurately.

With reflective surfaces or objects, the LED light can result in overexposed pictures or reflections.

- Turn the microscope camera or the object to be viewed until the picture is aligned as desired.

Make sure no objects touch the lens as they might damage it. Do not touch the optics (the lens will get dirty).

- The microscope camera can be fixed at a given position using the table stand included in the delivery; the locking bolt (3) fixes the position of the table stand.

Handling

Never use the product immediately after it has been taken from a cold room to a warm one. The condensation that forms might destroy your unit.

Allow the device to reach room temperature before switching it on. This may take several hours. The housing must not be opened! Any maintenance or repair work may only be carried out by authorised personnel.

The following unfavourable conditions are to be avoided when installing or transporting the product:

- Excessive humidity or damp
- Extreme cold or heat
- Dust or flammable gases, fumes or solvents
- Strong vibrations
- Strong magnetic fields such as those found in the vicinity of machinery or loudspeakers

When setting up/using the product, ensure that the cables are not bent, over-tightened or squeezed.

Cleaning

Check the lens of the microscope camera occasionally, clean it of dust and dirt. Use an appropriate lens cleaning system, e. g. a clean, soft brush. Do not press too hard on the optics/mechanics of the camera, since they may be damaged. Furthermore, you may cause scratch marks on the lens.

Disposal

Dispose of an unserviceable product in accordance with the relevant statutory regulations.

Technical Data

SensorCMOS, 1600 x 1200 pixels (2 MP)
Resolution image max. 2560 x 2048 pixels (only via software)
Resolution video max. 640 x 480 pixels
Image refresh rate 30 images per second
USB connectionUSB2.0
Operating voltage5V= (via USB)
Illumination4 white LEDs
Magnification90x to max. 500x at 20" (51 cm) monitor
Area of focusingup to 10mm
Dimensions (Ø x L)approx. 37 x 140 mm

These operating instructions are a publication by Conrad Electronic SE, Klaus-Conrad-Str. 1, D-92240 Hirschau (www.conrad.com).

All rights including translation reserved. Reproduction by any method, e.g. photocopy, microfilming, or the capture in electronic data processing systems require the prior written approval by the editor. Reprinting, also in part, is prohibited.

These operating instructions represent the technical status at the time of printing. Changes in technology and equipment reserved.

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© Copyright 2011 by Conrad Electronic SE.
V1_1111_01/AB